

Sulfo-SDTB

Product Description

cat.number : [UP35300A](#) , 50mg
Name: Sulfo-SDTB
Chemical Structure: Sulfosuccinimidyl-4-o-(4,4-dimethoxytrityl) butyrate
 (C₂H₅NO₂), Molecular weight: 665.60
Storage: +4°C (x)

Directions for Use

Protocole¹: Quantification of Amine Groups Using Sulfo-SDTB:

- A stock buffer consisting of 50mM sodium bicarbonate, pH 8.5 was prepared. Amine bearing surfaces were cut and weighed. Sulfo-SDTB (3mg) was weighed and dissolved in 1ml dimethyl formamide. After thorough mixing, the sulfo-SDTB solution was brought up to a total volume of 50ml with the stock sodium bicarbonate buffer (working sulfo-SDTB solution). Stock buffer (1ml) and 1ml working sulfo-SDTB solution were added to each tube and reacted for 40 minutes at room temperature on an orbital shaker. Segments were then removed and washed twice in 5ml of distilled water on an inversion mixer. Immediately following the wash, 2ml of a perchloric acid solution was added to each segment. Segments were reacted for 15 minutes on the inversion mixer. The reaction solution (1ml) was then removed and absorbance at 498nm was measured. Using the extinction coefficient for sulfo-SDTB and the segment weight, amine content (pmoles)/Ti segment (mg) was determined

Product Information

Sulfo-SDTB is used in order to determine amino groups on matrice⁽¹⁾. The reaction between sulfo-SDTB and amino groups results in stable bonds at alkaline conditions. The simple determination of amino group is due to 4,4'-dimethoxytrityl, (absorbance at 498nm; extinction coefficient at 70000M⁻¹cm⁻¹) which is released after the reaction of sulfo-SDTB on amino groups. The unbound reagent must be removed before acidic developing solution is added, therefore the applications of this assay are limited to testing solid phase supports. This assay procedure may be developed for amine estimation of soluble proteins by binding the protein of interest to microwell plates, which would allow the washing off of excess reagent without removing the protein.

Additional Information

Reagent for R&D *in vitro* use only

Literature

[Donald 2002](#): Development of a Bioactive Surface for Titanium Implant Devices Donald J. Dempsey, +Matthew D. Phaneuf, +Frank W. LoGerfo, +William C. Quist ; ©2002 SOCIETY FOR BIOMATERIALS 28th Annual Meeting Transactions
 Gaur 1989: Gaur, R.K and Gupta, K.C. (1989). A spectrophotometric method for the estimation of amino groups on polymer supports. Anal.Biochem.180, 253-258.

Contact uptima – Interchim for any question

Rev. C06E

Contact your local distributor

uptima@interchim.com

Uptima, powered by

 213 Avenue J.F. Kennedy - BP 1140
 03103 Montluçon Cedex - France
 Tél. 04 70 03 88 56 - Fax 04 70 03 82 66